

EJB-H series

Electronics balances

User's guide

Contents

SECTION 1 INTRODUCTION

SECTION 2 SPECIFICATIONS

SECTION 3 INSTALLATION

SECTION 4 KEY DESCRIPTIONS

SECTION 5 OPERATION

- 5.1 Zeroing the display
- 5.2 Taring
- 5.3 Weighing a sample
- 5.4 Percent weighing
- 5.5 Parts counting
- 5.6 Accumulated

SECTION 6 BATTERY OPERATION

SECTION 7 RS-232 OUTPUT

SECTION 8 PARAMETERS

SECTION 9 APPENDIX

- 9.1 Error code
- 9.2 Weighing unit listing


SECTION 1 INTRODUCTION

The EJB-H series of electronics balances provides an accurate, fast and versatile series of general purpose balances with counting, % weighing functions.

There are 4 models in each series, with capacities up to 3000g. They all have stainless steel weighing platforms on an ABS base assembly.

All the keypads are light touch switches and the displays are large easy to read liquid crystal type displays (LCD). The LCD's are supplied with a backlight and have loading parts counting.

All units include automatic zero tracking, automatic tare, and an accumulation facility that allows the count to be stored and recalled as an accumulated total.

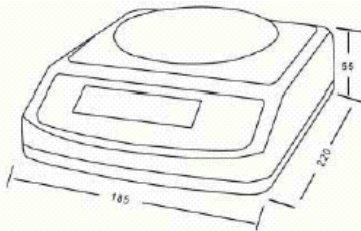

SECTION 2 SPECIFICATIONS

2.1 EJB-H series

EJB-H SERIES				
Model #	EJB-H-300	EJB-H-600	EJB-H-1500	EJB-H-3000
Max. Capacity	300 g	600 g	1500 g	3000 g
Readability	0.005 g	0.01 g	0.02 g	0.05 g
Resolution	1/60,000			
Tare range	100 g	300 g	600 g	1000 g
Min. Capacity	0.1 g	0.2 g	0.4 g	1 g
Linearity ±	0.01 g	0.02 g	0.04 g	0.1 g
Platter	Φ 120mm		136mm × 162mm	
wind shield	Yes			None
Units of Measure	g / ct / lb / oz / d / gn / ozt / dwt / mm / tl.T / tl.c / tl.t / t			

2.2 Common Specifications

Interface	RS-232 Output Optional
Stabilization Time	2 Seconds typical
Operating Temperature	0°C - 40°C / 32°F - 104°F
Power supply (external)	9V/800mA or built-in rechargeable battery
Calibration	Automatic External
ADC	Σ-Δ
Internal counts	500,000(EJB) or 1,000,000(EJB-H)
Display	20mm high 6 digits LCD with auto backlight and loading bar graph
Balance Housing	ABS Plastic, Stainless Steel Pan
Pan Size	Φ120mm / 136mm × 162mm
Overall Dimensions (wxdxh)	180mm x 220mm x 85mm
Gross Weight	2.6kg(include rechargeable battery)
Other Features and Specs	accuracy enhancement for parts counting and percentage weight function

SECTION 3 INSTALLATION

3.1 GENERAL INSTALLATION

The scales should be sited in a location that will not degrade the accuracy.

Avoid extremes of temperature. Do not place in direct sunlight or near air conditioning vents.

Avoid unsuitable tables. The tables or floor must be rigid and not vibrate. Do not place near vibrating machinery.

Avoid unstable power sources. Do not use near large users of electricity such as welding equipment or large motors.

Avoid high humidity that might cause condensation. Avoid direct contact with water.

Do not spray or immerse the scales in water.

Avoid air movement such as from fans or opening doors. Do not place near open windows.

Keep the scales clean.

Do not stack material on the scales when they are not in use.


3.2 INSTALLATION of EJB-H SERIES

Level the scale by adjusting the four feet. The scale should be adjusted such that the bubble in the spirit level is in the centre of the level and the scale is supported by all four feet. If the scale rocks readjust the feet.

Attach the adapter to the connector on the side of the scale. Please use the power as the label.

SECTION 4 KEY DESCRIPTIONS

Zero or

Set the zero point for all subsequent weighing. The display shows zero. A secondary function , of "Enter" key when setting parameters or other functions.

Tare or

Tares the scale. Stores the current weight in memory as a tare value, subtracts the tare value from the weight and shows the results. This is the net weight.

%

Enters the percent weighing function.

When in percentage mode and current value is zero, press this key to return normal weighing mode.

Print or

To print the results to a PC or printer using the optional RS-232 interface. It also adds the value to the accumulation memory if the accumulation function is not automatic.

Secondary function (**ESC**), is to return to normal operation when the scale is in a parameter setting mode.

MODE or

This key will select either kilograms, pounds, ounce for the weighing unit when weighing mode.

A secondary function, change current value for parameters or other functions. When counting mode, allows the weight, unit weight, and count to be seen when parts counting.

Smpl or

Use this key to enter counting mode when normal weighing mode, also use this key to return normal weighing mode when counting Mode.

ON/OFF or

Use this key to turn on or turn off the power.

SECTION 5 OPERATION

5.1 ZEROING THE DISPLAY

You can press the **ZERO/ENTER** key at any time to set the zero point from which all other weighing and counting is measured, within 10% of power up zero. This will usually only be necessary when the platform is empty. When the zero point is obtained the display will show the indicator for zero.

The scale has an automatic rezeroing function to account for minor drifting or accumulation of material on the platform. When the weight goes over 20 % of the full capacity, "read zero" function is obsolete.

5.2 TARING

Zero the scale by pressing the **ZERO/ENTER** key if necessary. The zero indicator will be on.

Place a container on the platform, a value for its weight will be displayed.

Press the **TARE** key to tare the scale. The weight that was displayed is stored as the tare value and that value is subtracted from the display, leaving zero on the display. The "TARE" indicator will be on. As product is added only the weight of the product will be shown. The scale could be tared a second time if another type of product was to be added to the first one. Again only the weight that is added after taring will be displayed.

When the container is removed a negative value will be shown. If the scale was tared just before removing the container this value is the gross weight of the container plus all product that was removed. The zero indicator will also be on because the platform is back to the same condition it was when the **ZERO/ENTER** key was last pressed.

5.3 WEIGHING A SAMPLE

To determine the weight of a sample first tare the empty container then place the sample in the container. The display will show the weight and the units of weight currently in use.

5.4 PERCENT WEIGHING

The scale will allow a sample weight to be shown as 100%. Then any other weight placed on the scale will be displayed as a percentage of the original sample. For example if 350g is placed on the scale and the % key is pressed the display will show 100.00%.

Removing the 350g weight and putting a 300g weight on the scale the display will show 85.71% as 300g is 85.71% of 350g.

Note: the scale may jump by large numbers unexpectedly if small weights are used to set the 100% level. For example if only 23.5g is on a scale with 0.5g increments and the scale is set to 100%, the display will show 100.00%, however a small change of weight will cause the display to jump to 102.13% as one scale division (0.5g) increase to 24.0g will be equivalent to a 2.13% increase.

Removing the 350g weight and putting a 300g weight on the scale the display will show 85.71% as 300g is 85.71% of 350g.

Pressing the % key will return the scale to weighing.

5.5 PARTS COUNTING

When the scale is showing weight, pressing the **SMPL** key will start the parts counting function.

Before beginning, tare the weight of any container that will be used, leaving the empty container on the scale. Place the number of samples on the scale. The number should match the options for parts counting, 10, 20, 50, 100 or 200 pieces.

Press the **SMPL** key to begin. The scale will show "SP 10" asking for a sample size of 10 parts. The display will cycle through the options: 10,20, 50, 100, 200 and back to 10.

Press the **SMPL** key again when the number matches the number of parts used for the sample. As more weight is added the display will show the number of parts (pcs).

Press the **MODE** key to display unit weight (g/pcs), Total weight (kg) or the count (pcs). Press the **SMPL** key to return to normal weighing.

5.6 ACCUMULATION

When the scale is set to manual accumulation the weight displayed will be stored in memory when the **PRINT** key is pressed and the weight is stable.

The display will show "ACC 1" and then the total in memory for 2 seconds before returning to normal. If the optional RS-232 interface is installed the weight will be output to a printer or PC.

Remove the weight, allowing the scale to return to zero and put a second weight on. Press **PRINT**, the display will show "ACC 2" and then show the new total.

Continue until all weights have been added.

To view the totals in memory press the **PRINT** key when the scale is at zero. The display will show the total number of items "ACC xx" and the total weight before returning to zero. The totals will also be printed via the RS-232 interface.

To erase the memory press **PRINT** to view the totals and then press the **SMPL** key to clear the memory.

SECTION 6 BATTERY OPERATION

When the battery needs change the indicator will turn on. Please change battery or use AC adapter after this indicator on.

The scales can be operated from the battery if desired. The battery life is approximately 50 hours.

To charge the battery simply plug into the mains power. The scale does not need to be turned on.

The battery should be charged for 12 hours for full capacity.

There is an LED to indicate the status of battery charging. When the scale is plugged into the mains power the internal battery will be charged. If the LED is green the battery has a full charge. If it is Red the battery needs charge and yellow indicates the battery is being charged.

As the battery is used it may fail to hold a full charge. If the battery life becomes unacceptable then contact your distributor.

SECTION 7 RS-232 OUTPUT

The EJB-H Series of scales can be ordered with an optional RS-232 output.

Specifications:

- RS-232 output of weighing data
- ASCII code
- 9600 Baud
- 8 data bits
- No Parity

Connector: 9 pin d-subminiature socket

- Pin 2 Output
- Pin 3 Input, not used at this time
- Pin 5 Signal Ground

EJB-H can connect with PC, remote display or mini printer, setup see section 8

SECTION 8 PARAMETERS

The scale has 8 parameters that can be set by the user plus a method of entering the calibration section.

To set parameters press the **MODE** key during self checking

The display will show the first function, "F1 UNT".

Pressing the **MODE** key will cycle through the other functions.

Pressing **ZERO** key will allow you to set the function. Use the **PRINT/ESC** key to leave a parameter unchanged.

Attached: When display shows "TECH", press ZERO key to enter, display shows "PIN", press SMPL, PRINT, TARE key, and press ZERO key to enter technical parameters setting mode.

Parameters table.

FUNCTION	DESCRIPTION
F1 UNT	Sets the which units want to use: g / ct / lb / oz / d / gn / ozt / dwt / mm / tl.T / tl.c / tl.t / t, use MODE to switch, use ZERO key to sure.
F2 EL	Set the backlight to be on, automatic or off, EL on: EL backlight always on EL Au: EL backlight automatically turn on EL off: El backlight always off Use MODE key to switch, use ZERO key to sure
F3 PRT	Set RS-232 P1 PRT connect with mini printer P2 RE connect with remote display or PC Use MODE key to shift, use ZERO key to sure
TECH	Pin Enter the programming and calibration menus by entering the correct password (SMPL, Print, TARE) and press ZERO key to enter.
P1 LIN	Liner calibrate After you press ZERO key, display will show PIN again, press PRINT, SMPL, % key and ZERO key to enter, display will show UNLOAD, remove all weight from the pan and then press the ZERO key to set zero point. The display will then show the first calibration weight request. Put this weight on the platform and then press the ZERO key when stable. The scale should be stable before pressing the ZERO key to accept a weight. The stability indicator will turn on to show the value is stable. After first weight calibrate, display will show the next calibrate weight, put calibrate weight like display shows, then press ZERO key. After all calibrate weight completed, press ZERO to complete the calibration. The scale will begin counting back to zero after the last weight has been selected.
P2 CAL	Normal calibrate After enter cal mode, the display will show "unLoAd". Remove all weight from the pan and then press the ZERO key to set the initial zero point. The display will then show the calibration weight request, you can use MODE key to select the calibrate weight. Put the weight you selected on the platform and then press the ZERO key when stable. The scale should be stable before pressing the ZERO key to accept a weight. The stability indicator will turn on to show the value is stable. The scale will begin counting back to zero after the last weight has been selected.
P3 CNT	This display will show xxxxxx for indicating the internal counts, you can press ZERO key to escape
P4 AZN	This option is used to select the auto zero maintain. Options : 0.5d, 1d, 2d, 4d, press ZERO key to sure, press PRINT key to escape

SECTION 9 APPENDIX

9.1 ERROR CODES

During the initial power-on testing or during operation it is possible the scale may show an error message. The meaning of the error messages is described below.

If an error message is shown repeat the procedure that caused the message, turning the balance on, calibration or other functions. If the error message still is shown then contact your dealer for further support.

ERROR CODE	DESCRIPTION	POSSIBLE CAUSES
Err 4	Initial Zero is greater than allowed (4% of maximum capacity) when power is turned on or when the ZERO/ENTER key is pressed	Weight on the pan when turning the scale on. Excessive weight on the pan when zeroing the scale. Improper calibration of the scale. Damaged load cell. Damaged Electronics.
Err 5	Keyboard Error.	Improper operation of the scale.
Err 6	A/D count is not correct when turning the scale on.	Platform not installed. Load cell damaged. Electronics damaged.

To view the A/D count press the **ZERO/ENTER** key and the % key at the same time while the scale is performing the initial check at power-on. Press **ZERO/ENTER** to return to normal.

9.2 WEIGHING UNIT LISTING

No.	Unit Mark	Unit Name	Conversion "g"
1	g	gram	
2	g	gram (hi resolution)	
3	ct	net carat	=0.2g
4	lb	lb	=453.59237g
5	oz	ounce	=28.349523125g
6	d	dram	=1.7718451g
7	gn	grain	=0.06479891g
8	ozt	troy ounce	=31.1034768g
9	dwt	Penny weight	=1.55517384g
10	mm	Momme	=3.749996g
11	tlT	tael tw	=37.49995g
12	Tlc	tael chn	=37.799375g
13	Tlt	troy tael	=37.4290018g
14	t	tola	=11.6638039g

“ENDEAVOUR” is an internationally reputed brand name in Electronic Weighing Instruments since last 11 years.

Our products are produced in India & Taiwan and supplied & serviced from our India, UAE and Kenya hubs; distributed through a world wide dealer network.

All “ENDEAVOUR” products are carrying advanced technology, designed by our R & D experts for specific commercial & non-commercial use in Laboratory, Medical, Business and Industrial environments.

Brief summary of “ENDEAVOUR” product range is...

- Laboratory Balances
- Jewelry Balances
- Pocket Scales
- Gem Scales (Carat Scales)
- Table Top Scales / Super Market Scales
 - Simple Weighing Scales
 - Price Computing Scales
 - Piece Counting Scales
 - Wash Proof/ Water Proof Scales
- Platform Scales
- Deck Scales
- Pallet scales
- Crane Scales
- Hanging (fishing) Scales
- Postal / Parcel Scales
- Bathroom Scales
- Kitchen Scales
- Baby Weighing Scales
- Truck Weighing Scales (Weighbridges)
- Weighing Indicators
- Load Cells
- Accessories
- Customized auto weighing systems

“ENDEAVOUR” operates on approved quality management system and is certified to ISO 9001: 2000

Visit our web : <http://www.dubaiweighing.com> for more details

CONTACT for any further assistance



PARAG TRADING CO., LLC

22, Bin Sabt Bldg., P-II Al musalla, P.O. Box # 51687, Bur-Dubai, DUBAI, U A E

Phone : +971-4-3522077 □ Fax : +971-4-3522088

Email : paragtrd@emirates.net.ae

© PARAG TRADING CO., LLC - 2004

©Copyright by PARAG TRADING CO., LLC All rights reserved.

No part of this publication may be reprinted or translated in any form or by any means without the prior permission of PARAG TRADING CO., LLC

PARAG TRADING CO., LLC reserves the right to make changes to the technology, features, specifications and design of the equipment without notice.

All information contained within this publication was to the best of our knowledge timely, complete and accurate when issued. However, we are not responsible for misimpressions which may result from the reading of this material.